



## Convert to csv file


1. To convert the spectrum to reflectance in csv file open the ASD program ViewSpecPro in the laptop of the ASD FieldSpec
2. Open the files to convert File>Open, select all files Ctrl+A and click 'Open'

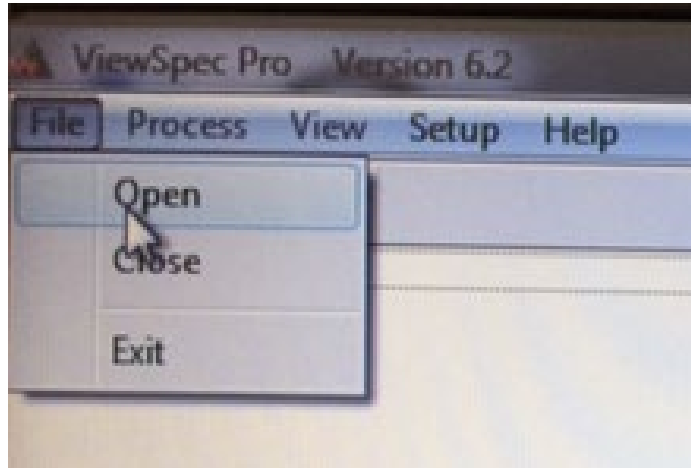

3. Select again all files that are in the screen, then go to Process>ASCII Export.

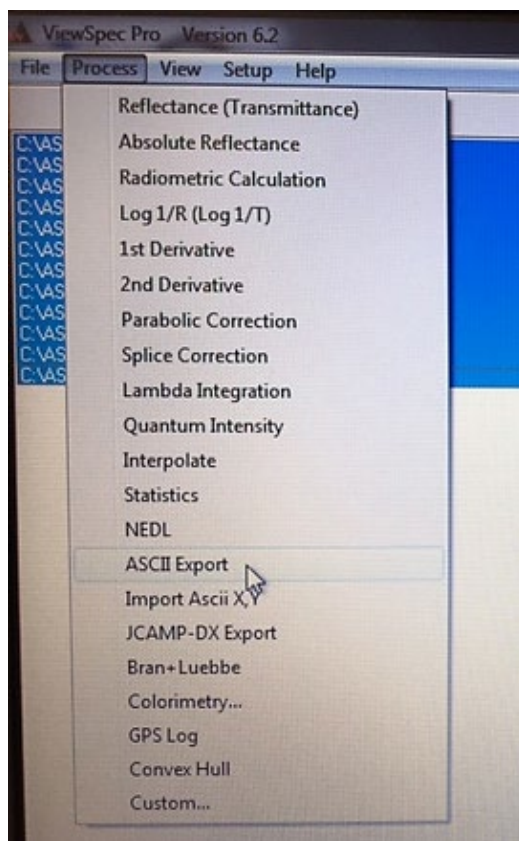

4. In the section 'Data Format for .asd files only' select reflectance. In the section 'Field Separator', select 'Comma' and tick 'output to a single file', click OK and save with 'filename.csv', select Open (See figure below). ViewSpecPro will convert the ASD files in one single csv file, which is a format that can be opened in excel.

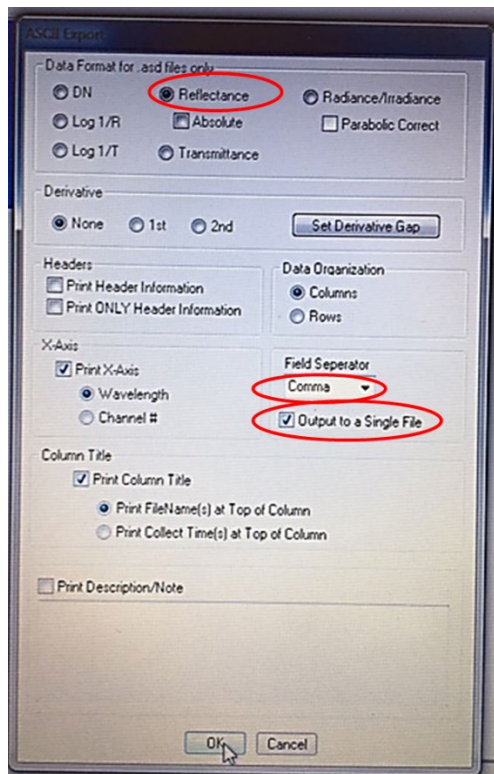

5. Copy the folder created at the beginning to a USB 'C:/ASD/Viri/ create a sub folder (YYMMDD\_your name)'.